


## Analysis and Data Management

The **Cytek® Amnis® IDEAS®** software is used to analyze data collected on the **ImageStream®** system. There are three types of image file: **Raw Image File (.rif)**, **Compensated Image File (.cif)**, and **Data Analysis File (.daf)**. FCS files can be analyzed in IDEAS® software or third party analysis software and will contain the feature data without the images.

### Opening files using the Start Analysis

1. Launch IDEAS® software.
2. Click **Start Analysis**, and browse for the appropriate **data file**. Or use “File” “Open” and choose a .rif, .cif, (skip to step 4) or .daf (skip to step 6). File extensions need to be visible in windows to know what the file type is.
3. Browse for an existing **compensation** matrix (.ctm file) or create a new one by clicking **New Matrix** and follow the steps outlined in the compensation QS guide.
4. Select an analysis **template** (.ast file) or a .daf with the analysis you wish to apply. If no template is chosen the default template will be used.
5. Select the active channels used in the experiment, and the **Image Gallery Display** properties will be applied to the images. They can be manually adjusted following the reporting QS guide.
6. Select the appropriate **wizard** for your application and follow the steps to completion. Basic Immunophenotyping can be done using the “**Begin Analysis**” wizard. Use the “**feature finder**” wizard if your application is not in the default list.
7. To add additional analysis use the **histogram** and **dot plot** icons in the tool bar, or select **Guided analysis** and choose a building block.
8. Add additional **Masks** as needed. (Analysis / Masks / New / Function).
9. Add additional **Features** as needed. (Analysis / Features / New).
10. When the analysis is complete **save** the data analysis file (.daf) and analysis temple (.ast) using the options in the file drop down list. This analysis can be applied to any data file following the steps above, or choosing “file” then “open” and applying the .ast and .ctm files.
11. When the wizard is complete, validate the regions by opening the control files using the saved analysis template and compensation matrix. Also verify the data in the statistics report under “reports” and “**define statistics report**”.
12. To **batch process** additional files, select tools/ batch data files and refer to the Batching Multiple Files QS guide 3.
13. To **report** the data, right click plots or images, and chose copy to clipboard. You can then paste the data into any third party reporting software.

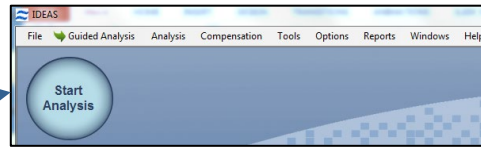

# IDEAS® Software Quick Start Guide

## To save an analysis template file:

1. From the File drop down option, Select **Save As Template File (.ast)**.
2. Enter the **name** of the analysis template to save.
3. Click **Save**. Any file can now be opened and analyzed using this template.

## To batch process data files:

1. From the **Tools** menu, select **Batch data files**.
2. Click **Add Batch**.
3. Click **Add Files**, and ctrl select all the data files in the experiment.
4. Select the **compensation** and **analysis** template to use.
5. Click OK, and when all critical batches have been added, select **Submit Batch**, to start the batch.

## To merge Raw Image Files:

1. From the **Tools** menu, select **Merge .rif Files**. The Merge .rif Files screen appears.
2. Click **Add Files**, and Ctrl-select the .rif files to merge.
3. Click **OK** and name the file with a unique file name.
4. Click **Save**.


## To create a new data file from populations:

1. From the **Tools** menu, select **Create Data File from Population**.
2. Check the **Raw Image File (.rif)** box to create a .rif file and click the browse button to select a name. Check the **Compensated Image File (.cif)** box to create a .cif.
3. Check the desired populations in the **Select population** list, and hit OK.

## To Export Image and Feature Data for multiple files:

1. Under the **Tools** menu, select the type and format of the data to export.
2. To export **Feature data**, select export features, for images select export .tif Images.
3. Select the population and features to export.
4. Select **Add files**, and Ctrl-select the data files to export.
5. Chose the format and population of data to export (i.e. FCS), and press OK.

## File Structure Overview (.rif, .cif, .daf)

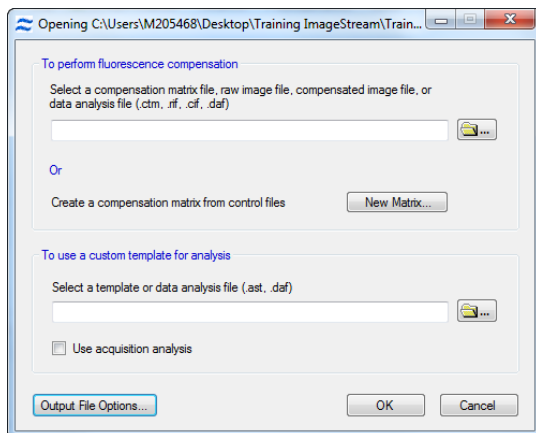


2.) Open a raw data file .rif

3.) Navigate to a compensation matrix.

4.) Navigate to an analysis template.

12.) Once analysis is validated **Batch** process experiment, and review statistics report.

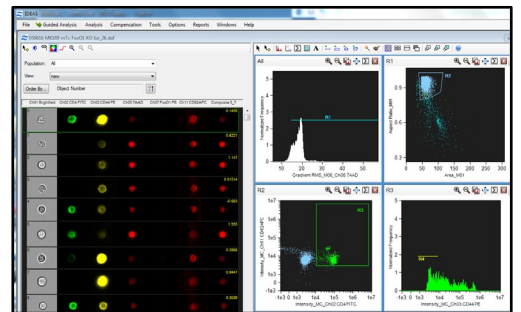

©2023 Cytek Biosciences, Inc. All rights reserved. Cytek, Amnis, IDEAS and ImageStream are trademarks of Cytek Biosciences, Inc. All other trademarks are the property of their respective owners.